

DLA - Proficiency Tests GmbH

Hauptstr. 80 23845 Oering/Germany

Tel: +49-(0)4532/9183358 Mob: +49-(0)171/1954375 Fax: +49-(0)4102/9944976 eMail: pt@dla-lvu.de Internet: www.dla-lvu.de

DLA, Hauptstr. 80, 23845 Oering/Germany

Datum / Date: 8. Mar. 2025

Dear participants,

Please find enclosed the material for the proficiency test (PT):

DLA ptALS1 (2025) - Allergen-Screening I - 4 Samples qualitative: Almond, Brazil Nuts, Cashew, Coconut, Hazelnut, Macadamia, Pecan, Pine Nuts, Pistachio and Walnut (NEW: gravimetric 2/3 and 1/6 values of new "Action Levels" for all allergens included)

There are 4 different samples possibly containing the allergenic parameters for qualitative determination. The allergens are spiked in a simple food matrix (rice flour / maltodextrin) and all parameters are present in 2 of 4 samples providing 2/3 and 1/6 of the new Action Levels (VITAL 4.0, 2023), respectively. The "action level" in mg/kg is calculated from the allergen protein reference dose per 100g food portion.

You are registered for

- a) 1 analysis method: or
- b) 2 or more analysis methods:)

Please note the attached information on the proficiency test.

<u>New:</u> Please enter your final results online in our <u>PT customer portal</u> **my DLA | participant's portal**. You will receive further information on this by e-mail, in particular about access to the portal.

Last deadline is April 2025-04-25.

After the deadline no results can be accepted.

We are looking forward to any suggestions or questions! We wish you a successful performance of the proficiency test!

Kind regards,

Matthias Besler-Scharf & Alexandra Scharf

On behalf of the DLA-Team

DLA - Proficiency Tests GmbH



Hauptstr. 80 23845 Oering/Germany

Tel: +49-(0)4532/9183358 Mob: +49-(0)171/1954375 Fax: +49-(0)4102/9944976 eMail: pt@dla-lvu.de Internet: www.dla-lvu.de

Information on the Proficiency Test (PT)

PT number	DLA ptALS1 (2025)
PT name	Allergen-Screening I - 4 Samples qualitative: Almond, Brazil Nuts, Cashew, Coconut, Hazelnut, Macadamia, Pecan, Pine Nuts, Pistachio, Walnut in simple food matrix
Sample matrix	Samples 1-4: Food matrix / ingredients: rice flour (appr. 75%), maltodextrin (appr. 25%), other food additives and allergenic foods
Number of samples and sample amount	4 different Samples 1-4: 20 g each
Storage	Samples 1 - 4: room temperature (PT period), cooled 2 - 10°C (long term)
Intentional use	Laboratory use only (quality control samples)
Parameter	Qualitative results - Total food items are spiked with following 2/3 and 1/6 of the Action levels (calculated as total protein): Almond, Brazil Nuts, Cashew, Coconut, Macadamia, Pecan, Pine Nuts, Pistachio, Walnut (as Total Protein 6,7 mg/kg and 1,7 mg/kg, each) Hazelnut (as Total Protein 20 mg/kg and 5,0 mg/kg)
Methods of analysis	The analytical methods ELISA (+ Lateral Flow), PCR, NGS and LC/MS can be applied for qualitative determinations.
Notes to analysis	The analysis of PT samples should be performed like a routine laboratory analysis. In general we recommend to homogenize a representative sample amount before analysis according to good laboratory practice, especially in case of low sample weights.
Result table	One result each should be determined for Samples 1-4. The results should be filled in the result entry table.
Units	positive / negative (limit of detection mg/kg)
Number of digits	at least 2
Result submission	online via my DLA participant's portal (https://my.dla-pt.com) you will receive further information about the access by e-mail
Last Deadline	the latest <u>April 2025-04-25</u>
Evaluation report	The evaluation report is expected to be completed 6 weeks after dead- line of result submission and will be provided as a PDF file in the DLA Participant Portal (https://my.dla-pt.com/).
Coordinator and contact person of PT	Matthias Besler-Scharf PhD

^{*} Control of mixture homogeneity and qualitative testings are carried out by DLA. Any testing of the content, homogeneity and stability of PT parameters is subcontracted by DLA.

page 2 of 2